

**Protect hygiene where
it matters most.**

kikkoman
Kikkoman Biochemifa Company



Easy Plate™

Prepared media
for microbiological
testing



MICROVAL®  nin

Pros

- Simple operation, results
- Reliable
- AOAC PTM, NordVal & MicroVal Certificate
- Ready to use, inoculate the sample and incubate.
- No need for special tools
- Reduced waste, incubation space and time
- Brightly colored colognes that are easily distinguishable
- High accuracy and correlation
- Fast Counting and Differentiation Software



HyServe

HyServe GmbH
Ammerthalstraße 7
Tel. +49 (0) 89 3703 1223
info@hyserve.com


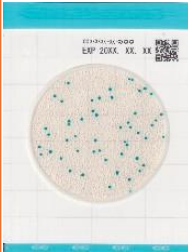
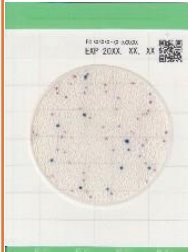



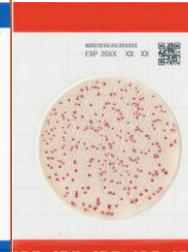
D- 85551
Kirchheim/München
Fax +49 (0) 89 3703 1225
www.hyserve.com

Protecting hygiene where it matters most

kikkoman
Kikkoman Biochemifa Company

A solution that facilitates microbiological testing!

Save time, labor, and money with Easy Plate. Easy Plate is a convenient, pre-prepared film media for faster and easier microbiological testing. All Easy Plates are AOAC validated and come in an easy-to-use, stackable film plate format that saves space, reduces waste, eliminates setup time and labor, and streamlines your testing procedures. Each Easy Plate is specialized for a specific target organism (or group) and includes options for aerobic bacteria (AC), coliforms (CC), E. coli/coliforms (EC), and Staphylococcus aureus (SA). More parameters will be available soon.

AC	CC	EC	YM-R	SA	EB	AC-R
Aerobic bacteria	Coliform	E. Coli / Coliform	Yeast and Mold Rapid	Staphylococcus aureus	Enterbacteriaceae	Aerobic Bacteria Rapid
						
35± 1°C			25± 1°C	35± 1°C / 37± 1°C		35± 1°C
48± 2 hours	24± 1 hours	24± 1 hours	48± 2 hours	24± 1 hours	24± 1 hours	24± 1 hours
Storage (Unopened): 2-8°C						
18 Months after danufacturing date				12 Months after manufacturing date.		
AOAC RI PTM Certified / MicroVal Certified / Nordval Certified					In process	
Packing: 100 tests (25 tests x4 Bags)						

Shelf life and storage

Storage conditions: 2-8°C (do not freeze)


4°C: 3 months (unopened)

25°C: 14 days (unopened) 30°C: 5 days (unopened)

After opening the package, fold the open end of the package at least 2 times and tape it down.

Easy to use


Due to the unique hydrophobic ring around the plate, it is ensured that the sample suspension is evenly distributed over the entire surface of the plate without the need for a spreader.

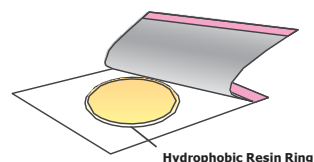


+ Dilution rack
PBS / MRD

+ ProMedia ST-25 PBS
Sampling Swab

Save Space

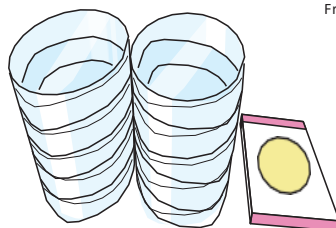

Amount of CO₂ Emitted

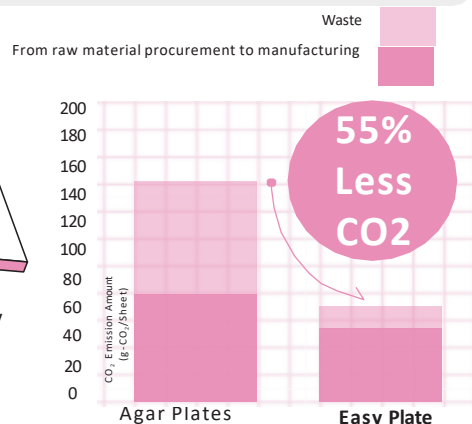

HyServe

HyServe GmbH
Ammerthalstraße 7
Tel. +49 (0) 89 3703 1223
info@hyserve.com

D- 85551
Kirchheim/München
Fax +49 (0) 89 3703 1225
www.hyserve.com