

**Fast and Effective
Allergen Testing**

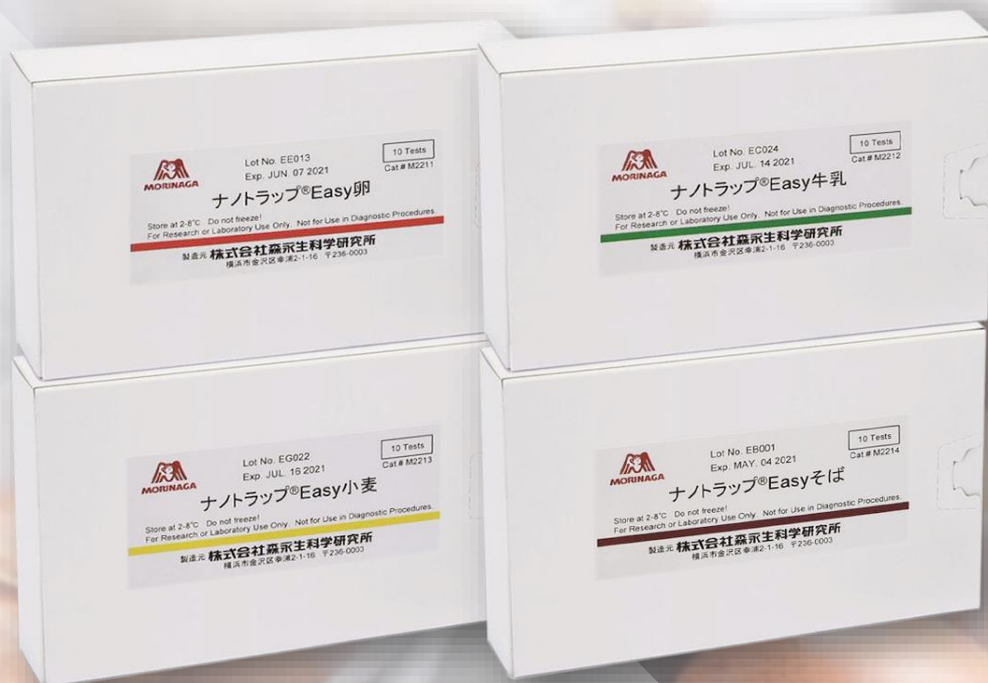


RAPID TEST

For food allergens

EASY

Make it
your
product:
OEM



Characteristics


- Simple and easy to use
- Specialized for rinse and swab water
- No sample preparation time required
- No special equipment needed
- Fast results in 10 minutes


HyServe


HyServe GmbH
Ammerthalstraße 7
Tel. +49 (0) 89 3703 1223
info@hyserve.com

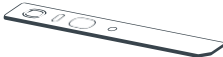
D- 85551
Kirchheim/München
Fax +49 (0) 89 3703 1225
www.hyserve.com

For Swab Samples

- 

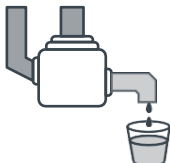
Sample the specific area
- 


Shake vigorously
- 

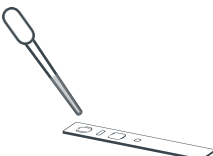
Add 4 drops to the test strip
- 


Incubate for 10 minutes at room temperature.

For Water Sampling

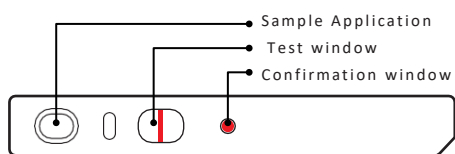
- 

Collect the water sample
- 

Designate it as a test solution
- 

Add 4 drops to the test strip
- 

Incubate for 10 minutes at room temperature.



Positive: A red violet line in a test window along with the red color in a confirmation window represents a positive allergen test.

Item-No.	Kit Specifications	Measurement Time	Limit of Detection (μg protein/ml)
1500 012	Egg (Ovomucoid)	10 minutes	0.5 ppm
1500 013	Casein (Casein)	10 minutes	0.5 ppm
1500 014	Gluten (Gliadin)	10 minutes	0.4 ppm
2 500 017	Soya (β -conglycinin)	10 minutes	0.5 ppm
2 500 018	Peanut (Multiple proteins)	10 minutes	0.5 ppm
2 500 019	Crustaceans (Tropomyosin)	10 minutes	0.5 ppm
2 500 020	Buckwheat (Multiple Proteins)	10 minutes	0.5 ppm

Conservation and shelf life

Conservation:

2-30°C (35 to 86°F), do not freeze.

Expiry date: Scope of delivery:

24 months (after manufacture)

1 x Box with 10 tests

Kit Components: Test Strip: 10 packs, Swab Kit: 10 bottless, Insert: 1 Sheet

HyServe

HyServe GmbH
Ammerthalstraße 7
Tel. +49 (0) 89 3703 1223
info@hyserve.com

D- 85551
Kirchheim/München
Fax +49 (0) 89 3703 1225
www.hyserve.com