




FOOD ALLERGEN

Rapid Test Easy

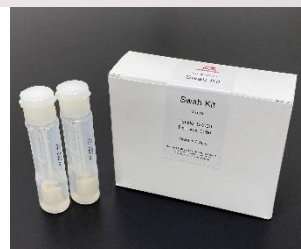

Kits

- Egg (Cat. # M2241)
- Casein (Cat. # M2242)
- Gluten (Cat. # M2243)
- Buckwheat (Cat. # M2244)
- Peanut (Cat. # M2245)
- Soya (Cat. # M2246)
- Crustacean (Cat. # M2247)
- Sesame (Cat. # M2248)
- Mustard (Cat. # M2249)

Features

- Simple and easy to use
- Suitable for surface swabbing and rinse water
 - * With the Easy for Sesame (M2248), only surface swab testing is available.
- No need for special equipment (swab kit included)
- 24-month shelf life
- Room temperature storage (2-30°C)

Included Kit Components

- Test Stick (10 packs)
- Swab Kit (10 bottles)
- Instruction

Features

Performance

Sample preparation time	Not needed
Assay time	10 min
Limit of detection for swab	0.5 ppm (μg protein / mL)
Limit of detection for rinse water	0.5 ppm (μg protein / mL)



Morinaga BioScience, Inc.
2-1-1 Shimosueyoshi,
Tsurumi-ku, Yokohama-shi,
230-8504
Japan



<https://www.morinaga-biosci.com>



info_miobs_e@morinaga.co.jp




TEL +81-45-586-2515
FAX +81-45-586-2517

SAMPLE COLLECTION


Surface Swabs



Swab the area of interest

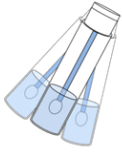


Shake vigorously

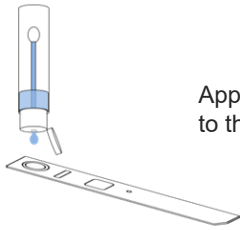


Apply 200 μ L (4 drops) to the test stick



Rinse Water

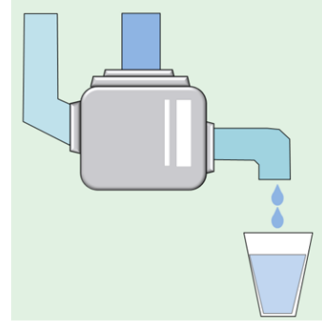

Collect rinse water

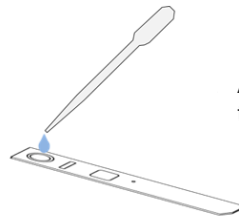

Apply 200 μ L (4 drops) to the test stick



Wait 10 min

RESULT



Negative

Positive

Explanatory Videos

Videos available in our website and YouTube channel show all the kit components and step-by-step sample preparation and assay procedure.

Related Food Allergen Products

QC Materials (Egg, Milk, Wheat)
Rapid Test Pro II
ELISA Kit II

Morinaga BioScience, Inc.

<https://www.morinaga-biosci.com>