Compact Dry X-SA medium for *Staphylococcus aureus*

40 plates/Platten/plaques/placas/lastre/placas

ID-No. 1 002 960

240 plates/Platten/plaques/placas/lastre/placas

ID-No. 1 002 961

500 plates/Platten/plaques/placas/lastre/placas

ID-No. 1 402 961

1400 plates/Platten/plaques/placas/lastre/placas

ID-No. 1 502 990

English

Deutsch

Français

Compact Dry X-SA is a ready to use, chromogenic plate for detection of Staphylococcus aureus

Sample pretreatment

Viable count in water or liquid foodstuff
Drop 1 ml of specimen (dilute if necessary) on the middle of the Compact Dry plate.

Viable count in solid foodstuff

Add buffer solution to the sample and homogenize by stomacher®. Drop 1 ml of specimen (dilute if necessary) in the middle of the dry sheet of the Compact Dry plate.

Viable count in swab test specimen

Use the swab to wipe the surface, put into the device with wiping solution. Drop 1 ml of wiping solution (dilute if necessary) in the middle of the Compact Dry plate. It is recommended to use "Swab for Compact Dry" offered by HyServe Id-No. 1 002 952/3 / 1 402 954 (40/240/600 pieces).

Test instructions

1. Open the cap and drop 1 ml of specimen in the middle of the Compact Dry plate.
2. Specimen diffuses automatically and evenly into the sheet and transforms the dried sheet into a gel within seconds.
3. Put the cap again on the plate and write the information needed on the memorandum section.
4. Turn over the capped plate and put in the incubator.
5. After incubation count the number of blue colonies. White paper placed under the plate helps to count the colonies.

Incubation time 24 hours ± 2
Incubation temperature 35 ± 2 °C
AOAC approved 35 ± 2 °C
NordVal/MicroVal approved 37 ± 1 °C

Interpretation of the results

Blue Colonies are *Staphylococcus aureus*.

Storage and shelf life

Keep at room temperature (+ 5 to +30 °C). Total shelf life 21 months after manufacturing.

Notes

- High cell concentrations on plates will cause the entire growth area to become colored/bluish. In this case dilute the sample.
- After use please follow the current disposal regulations.
- The growth area is 20 cm². The back of the plate has a grid carved of 1 cm x 1 cm to make the colony counting easier. In case of any difficulties to count colonies due to large number of colonies grown, total viable count can be obtained by multiplying 20 by an average number of colonies per grid counted from several grids.
- Compact Dry plates are produced at an ISO 9001 certified site.

- AOAC Approval No. 081001
- MicroVal approval No. 2008-LR14
- NordVal Certificate No. 042, ISO 6888-1:1999

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